


3D Biotek Cell Seeding Flow-Chart

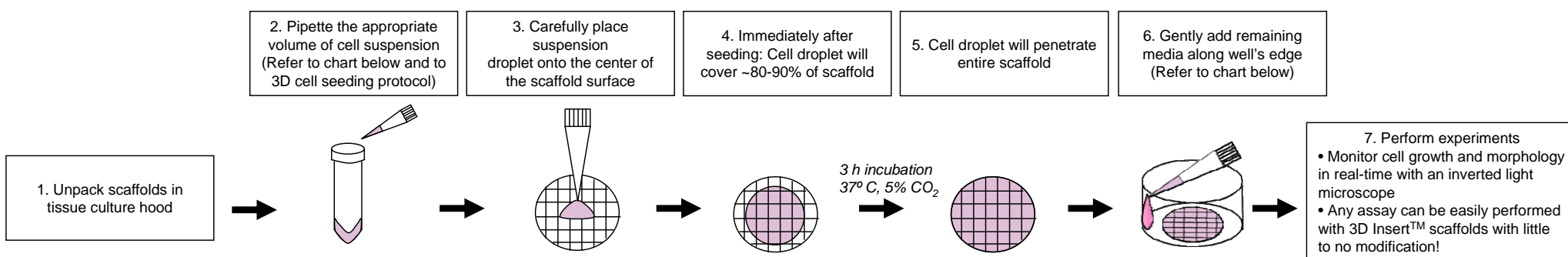

Scaffold Size	Seeding Volume (μl)			
	PS1520	PS3040	PCL3030	PCL3050
96-well	15	20	20	25
48-well	25	50	50	70
24-well	60	120	130	160
12-well	120	240	270	340
6-well	300	600	690	880

WARNING:
DO NOT touch scaffold with pipette tip!
DO NOT let droplet touch sides of wells.
This will decrease seeding efficiency!



Scaffold Size	Volume to Add After 3h (μl)			
	PS1520	PS3040	PCL3030	PCL3050
96-well	185	180	180	175
48-well	225	200	200	180
24-well	440	380	370	340
12-well	1380	1260	1230	1160
6-well	1700	1400	1310	1120

For further support, please visit us at <http://www.3DBiotek.com>

Technology Center of New Jersey ■ 675 US Highway One, North Brunswick, North Jersey 08902, USA ■ E-mail: info@3dbiotek.com